DNA Ladders

Omega Biotek offers a wide range of products for the accurate size and mass determination of DNA. Most DNA Ladder includes one or two higher-intensity bands for easy identification and orientation. The ladders are supplied premixed with loading buffer and are stable at room temperature for a few weeks. There is no need to heat, mix, or dilute ladders prior to loading them on gel.

100bp DNA Ladder plus

- CAT.No.: M08-01
- Packing: 250ul
- Concentration: 0.136 ug/ul
- Store: at 4° C for periods up to 6 months. For longer periods store at -20° C.
- Description: 100 DNA Ladder plus is designed for sizing and approximate quantification of wide range linear double-stranded DNA fragments on agarose gel. The ladder is composed of thirthy purified individual DNA fragments (in base pairs):3000,2000,1500,1200,1000,900,800,700, 600,500,400,300,200,100. It contains two high-intensity bands (1200bp and 500 bp) for easy identification and orientation.
- Direction:
 - \diamond This product can load directly, since it contains 1×loading buffer. It is free of fluorescence dye such as EtBr.
 - For a loading well less than 6mm in width, 5ul of this product is enough for clear bands. For a wider well, more product may be required.
 - ♦ Gel and buffer may effect the electrophoresis. For better result, please use high quality agarose and fresh running buffer. 2% agarose gel is recommended.

• Electrophoresis Map:

